

MSAK827

USB Powered LED Backlight


User Guide

Operation:

- Plug one end of the included micro usb cable into the USB Power Input and plug the other end into a 5V DC USB power supply (not included) or into the USB port of your computer.
- The backlight will automatically turn on when connected to a power source.
- For best results, turn off the built-in lights on your Dino-Lite digital microscope while using the backlight.

Safety Tips:

- Keep the backlight away from extremely hot or cold environments.
- Do not submerge the backlight in water.
- Avoid dropping the backlight as small pieces can break off and pose a choking hazard for pets and small children.
- If the backlight or USB cable becomes damaged, discontinue use immediately to avoid risk of electric shock.

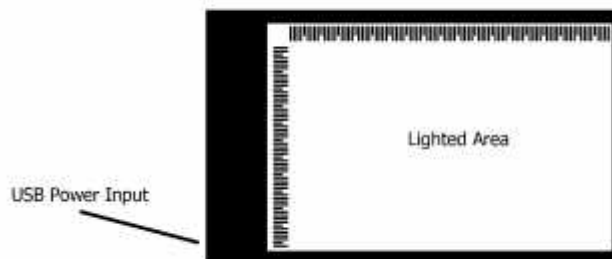

Specifications:

Frame size (mm): L148 X W105 X H5

Lighting area (mm): L116 X W92

Lighting Panel: Acrylic

Light Source: LED

Voltage: DV 5V

Rated Power: 3W

Power Port: Micro USB

Net Weight: 0.1kg